

AccuPower® RT-PCR PreMix

ISO 9001 Certified

AccuPower® RT-PCR PreMix is covered by Korea Pat. No. 162770 and US Pat. No. 5861251 & 6153412. Headquarters 49-3, Munpyeong-dong, Daedeok-gu, Daejeon 306-220, Korea

Fax:+82-42-930-8600

Seoul Office 1 floor, Donggwan Building 1566-2, Seocho-dong, Seocho-gu, Seoul 137-874 Korea Phone:+82-42-936-8500 Phone:+82-2-598-1094 Fax:+82-2-598-1096

Bioneer Inc. 1000 Atlantic Avenue Alameda, CA 94501 USA Toll free : 1-877-264-4300 Fax: 1-510-865-0350 E-mail:order.usa@bioneer.com

Order Korea Market only:1588-9788 E-mail: order@bioneer.co.kr Website: www.bioneer.com

AccuPower® RT/PCR PreMix contains all the components necessary for cDNA synthesis and amplification in one tube.

Advantages

Speed Substantial reduction in reaction setup time.

Stability As each tube of AccuPower® RT/PCR PreMix contains a stabilizer (patented in US and Korea), which can maintain the stability of the

RTase and DNA polymerase up to 2 years at -20°C. Pre-denaturation during PCR has no effects on the activity of the DNA

polymerase.

Reproducibility and Yield

The strict functional QC assays demonstrated highly consistent and reproducible RT/PCR performance. In most applications an

increase in yield is observed as compared to the standard reactions.

Simplicity The fewer manual steps allow reduction in potential errors. Each tube contains tracking dye and precipitant

for agarose gel electrophoresis, eliminating the needs for a separate loading buffer.

Experimental Protocol

· Mix the template RNA and the reverse primer in a sterile tube as indicated below;

Concentration of template RNA and primer

	Reaction volume	20 μl reaction	50 μl reaction		
Template	Total RNA	0.5~1.0 μg	1.0~2.0 μg		
RNA	Poly(A) RNA	0.05~0.1 μg	0.1~0.2 μg		
Primer	Primer	10~30 pmole	20~50 pmole		

- Incubate the mixture at 70°C for 5min and place it on ice.
- Transfer the incubated mixture and the forward primer to AccuPower® RT/PCR PreMix tube, then fill up the reaction volume with DEPC-DW.
- Dissolve the lyophilized blue pellet by vortexing, and briefly spin down.
- · Add mineral oil to each tube (This step is unnecessary when using a thermal cycler with top heating.).
- · Perform cDNA synthesis reaction as follows:

42°C, 60 min. (cDNA synthesis)

94°C, 5 min. (RTase inactivation)

• Perform PCR cycles according to the PCR condition.

(Annealing temperature and time need to be optimized for each primer/template combination.)

Trademark

AccuPower is a trademark of Bioneer Corporation.

Ordering Information

Tube type	Reaction	Cat. No	Description	Tube type	Reaction	Cat. No	Description
0.2ml Tube	20 ul	K-2055	0.2 ml thin-wall 8-strip tubes with attached cap / 96 tubes	96-well	20 ul	K-2262-4	thin-wall 96-well flat plate
		K-2055-B	0.2 ml thin-wall 8-strip tubes with attached cap / 480 tubes			K-2262-5	thin-wall 96-well full-skirted plate
	50 ul	K-2057	0.2 ml thin-wall 8-strip tubes with attached cap / 96 tubes			K-2262-6	thin-wall 96-well semi-skirted plate
		K-2057-B	0.2 ml thin-wall 8-strip tubes with attached cap / 480 tubes		5 ul	K-2084-1	thin-wall 384-well full-skirted plate
96-well	10 ul	K-2262-1	thin-wall 96-well flat plate	384-well	10 ul	K-2084-2	thin-wall 384-well full-skirted plate
		K-2262-2	thin-wall 96-well full-skirted plate		20 ul	K-2084-3	thin-wall 384-well full-skirted plate
		K-2262-3	thin-wall 96-well semi-skirted plate	0.5ml Tube	50ul	k-2056	0.5 ml thin-wall tubes with attached cap / 100 tubes

A complete product list appears on our web site at www.bioneer.com